


Selection and screening of stable mammalian cell lines

Stable cell lines are essential tools for genetic and phenotypic studies. To establish such lines, the transgenic cells which carry a defined genotype or exhibit a desired phenotype must be enriched or isolated from the parental population using a selectable marker. These markers typically encode either survival advantages such as an antibiotic resistance or detectable traits such as a fluorescence marker.

Selection can occur in a single step by applying selective pressure through antibiotics, which kill non-transgenic cells (*Alternative A*). Alternatively, fluorescence or surface markers allow rapid enrichment through screening and physical isolation (*Alternative B*). Refer to [SOP0036](#) for banking and tracking of cell lines.

This is a bench card. Full protocol available online.

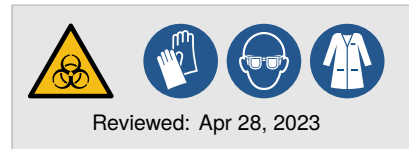

Procedures

A > Antibiotic selection of mammalian cell lines

Culture medium with antibiotics

- (1.) Determine the minimum effective antibiotic concentration for selection using a matrix titration format.
 - Prepare a 96-well plate with 100 μL of culture medium in all wells.
 - Perform a row-wise dilution of the antibiotic: Add 100 μL of the highest concentration to row A, and perform 1:2 serial dilutions down to row G. Leave row H as a no-antibiotic control.
 - In each row, seed column 1 with 100 μL of a $4.0 \times 10^4 \text{ mL}^{-1}$ cell suspension. Mix gently and transfer 100 μL to column 2. Continue serial dilution across columns to column 12. Discard excess.
 - Incubate the plate under standard conditions.
 - Assess cell survival under the microscope after 3–7 d.

- (2.) Select cells using the lowest antibiotic concentration that resulted in complete killing of non-transgenic cells. Include a negative control (non-transfected) to verify selection stringency. Useful ranges:

Cell line	Blasticidin S	G418 (Geneticin)	Hygromycin B	Puromycin	Zeocin
Typical duration	10 days	14 days	4 days	4 days	10 days
293T	2–10 $\mu\text{g}/\text{mL}$	Not determined	Not determined	1.0–5.0 $\mu\text{g}/\text{mL}$	Not determined
HeLa	5–10 $\mu\text{g}/\text{mL}$	Not determined	Not determined	1.0–2.0 $\mu\text{g}/\text{mL}$	Not determined
MCF 10A	2–10 $\mu\text{g}/\text{mL}$	150–400 $\mu\text{g}/\text{mL}$	100 $\mu\text{g}/\text{mL}$	0.5–2.5 $\mu\text{g}/\text{mL}$	750 $\mu\text{g}/\text{mL}$
iPS	10–20 $\mu\text{g}/\text{mL}$	Not determined	Not determined	Not determined	Not determined

- (3.) Maintain selection by feeding cells regularly. Expand once resistant colonies are visible.
- (4.) *Optional:* Perform a second round of limiting dilution under selection to ensure the selected cell population is monoclonal.

B > Screening of mammalian cell lines

- (1.) Physically separate the population into clones using limiting dilution, pipette-based cell picking, or flow cytometry (FACS).
- (2.) Feed cells regularly and monitor for clonal outgrowth. Split as needed.

Critical: In 96-well plates, border wells lose volume faster than center wells. Top up edge wells regularly to prevent desiccation.

Selection and screening of stable mammalian cell lines

(3.) Analyze each clone for expression of the transgene or phenotypic marker.



 [Recipe \(available online\)](#)  [Resources \(available online\)](#)  [Troubleshooting \(available online\)](#)  [Notes \(available online\)](#)

Lab Protocols — © B. C. Buchmuller. Licensed CC-BY-SA-4.0. Research use; provided as-is, without warranty. Current when printed. Visit <https://benjbuch.github.io/check/?q=698664f> or scan the QR code to read the full protocol or to check for updates.

